STERILE R
Sterile fluid path
Sterile Flüssigkeitsbahn
Chemin pour liquides stérile

Do not resterilize.
Nicht erneut sterilisieren.
Ne pas restériliser.

Buffy coat container (for blood stem cell rich fraction or whole blood)
Buffy-Coat-Beutel (für blutstammzellreiche Fraktion oder Vollblut)
Poche de couche leucocytaire (pour fraction riche de cellules souches hématopoïétiques ou pur-sang)

freeze volume (blood stem cell rich fraction or whole blood)
Einfriervolumen Buffy Coat (blutstammzellreiche Fraktion oder Vollblut)
volume à congeler de couche leucocytaire (fraction riche de cellules souches hématopoïétiques ou pur-sang)

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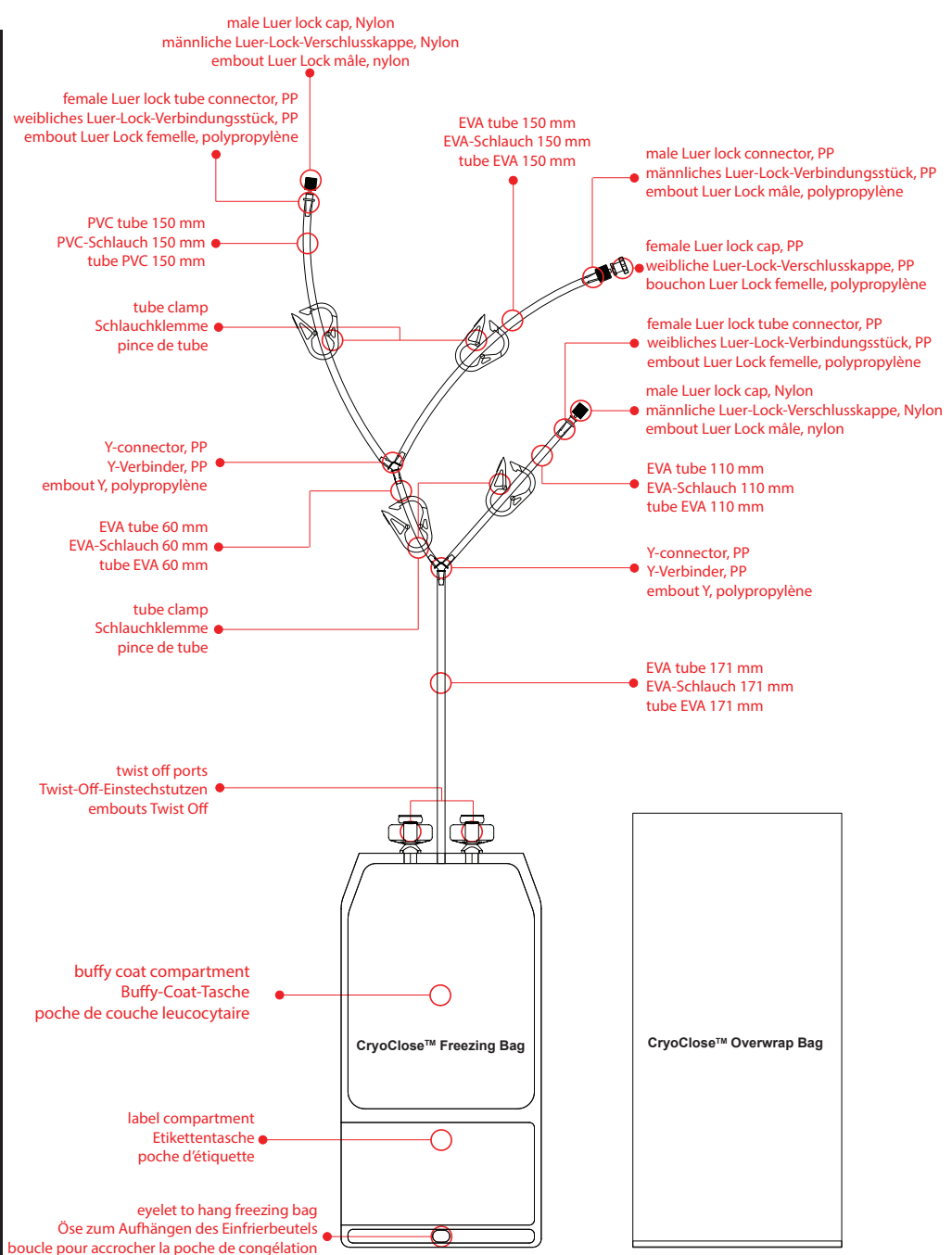
PHT DEHP

The yellowish coloured PVC-tube contains the phthalate DEHP (diethylhexyl phthalate). In order to prevent the release of DEHP into the blood which passes through the PVC tube respectively in order to reduce it to a minimum, anticoagulated blood shall be administered into the freezing blood bag via the PVC tube in less than one minute. For the same reason no DMSO, be it diluted or undiluted, shall pass through the PVC tubing. For the administration of DMSO into the freezing blood bag the transparent non-coloured EVA tubes shall be used. Repeated exposition to high doses of DEHP may exert nephrotoxicity or reproductive toxicity, especially in newborns and children or chronically ill patients such as for example dialysis patients or patients undergoing ECMO (extracorporeal membrane oxygenation).

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The maximum capacity is derived from the blood bag norm ISO 3826-1. It does not have any relevance for the intended use of the CryoClose™ Freezing Bag. For the intended use the 'recommended freeze volume' is pivotal only.

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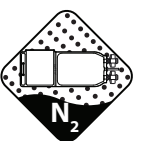
Intended use: The CryoClose™ Freezing Bag is intended to be used as a freezing container for cell suspensions, for cryopreservation at temperatures down to -196 °C in the vapour phase above liquid nitrogen. CryoClose™ Freezing Bag is intended for single use only. It is provided sterile within a sterile barrier peel bag. It is not to be re-sterilized. The fluid path of the CryoClose™ Freezing Bag is provided sterile and non-pyrogenic. Each CryoClose™ Freezing Bag is provided with a suitable overwrap bag.

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Preconditions for use of CryoClose™ Freezing Bag and general information

The CryoClose™ Freezing Bag is only intended to be used by specialized personnel, who have been trained in the cryopreservation of transplants of human hematopoietic stem cells and who dispose of the necessary experience. The processes of filling, closing, rate freezing, cryostorage as well as thawing when applying the CryoClose™ Freezing Bag are to be validated by the applying user.

Every box (transport packaging) that is labelled with a box label contains 28 CryoClose™ Freezing Bags. Each CryoClose™ Freezing Bag is accompanied by a suitable overwrap bag and sealed into a respective sterile barrier pouch. Each of these 28 sterile barrier peel pouches has been sealed into an outer transparent bag under clean room conditions, in order to facilitate the transport of the products into a cleanroom (see below). Caution: The transport box must not be opened with a knife to avoid damaging the enclosed packaging!

Transfer into clean room environment

28 CryoClose™ Freezing Bags are packed in a carton. These 28 CryoClose™ Freezing Bags have been individually packed into a transparent outside bag under clean room conditions. Before transfer into the clean room the transparent outside bag should be wiped clean and disinfected. After introduction into the clean room the transparent outside bag can be removed. The sterile barrier peel pouch should only be opened ultimately prior to application of the freezing blood bag, if possible under sterile conditions.

Before application the CryoClose™ Freezing Bag should be inspected visually: Do not use if package is damaged.

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